

# Mercury Collector

- **Collecting Mercury**
- **Safe storage**
- **Recycling and proper waste**


Mercury is toxic and has to be immediately collected after spilled in order to prevent health risks. Spilled mercury spreads immediately and is difficult with standard instruments. Using Mercury Collector from Ludwig Schneider it is extremely easy to collect and reuse spilled mercury. The mercury collector is composed of a Polyethylene.

Ludwig Schneider offer to reuse spilled Mercury from Thermometers or take care on a safe and environmental friendly waste of mercury.

Within the Mercury Collector screw cap lid there is a special plastic foam insert. On pressing this foam gently on the spilt mercury, the foam pores are opened and the mercury globules are being absorbed. On screwing the lid on the tube, the globules are released and dib through the perforated separation into lower part of the tub.

Mercury Collector from Ludwig Schneider will be supplied with instruction manual.

Mercury Collector from Ludwig Schneider Article No.: 61208

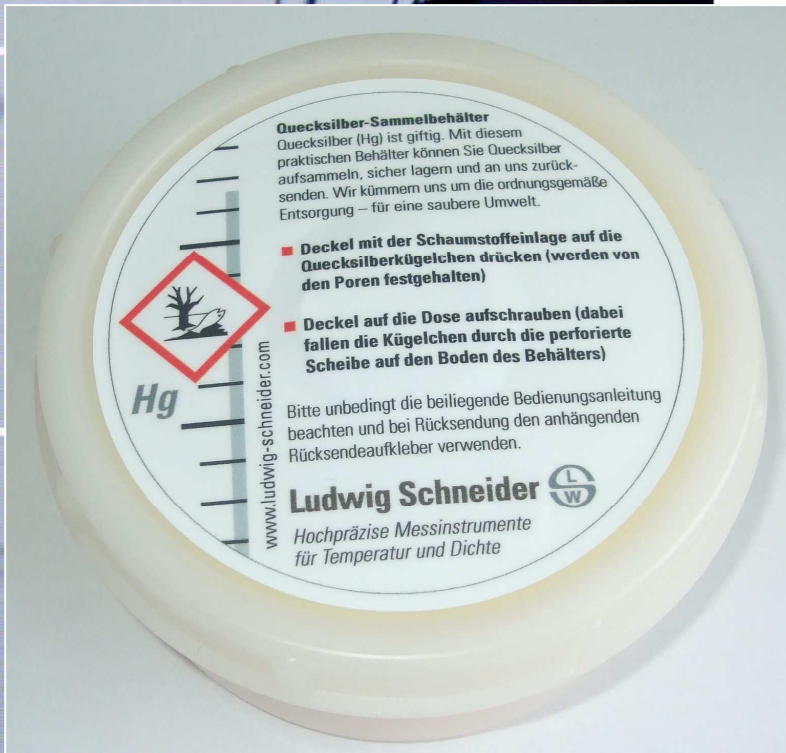

**Ludwig Schneider** 

**Ludwig Schneider GmbH & Co. KG**

Postfach 15 61 · 97865 Wertheim

Am Eichamt 4 · 97877 Wertheim

Tel.: +49-93 42- 85 60-0 · Fax: +49-93 42-8 46 71

e-Mail: [info@ludwig-schneider.de](mailto:info@ludwig-schneider.de)

[www.ludwig-schneider.com](http://www.ludwig-schneider.com)